

## Moltox EC Tri PC Plates (21-199)



**Intended Use:** 21-199 plates are used for the phenotypic confirmation of *E. coli* WP2 strains; tests for tryptophan auxotrophy and pKM101 presence (pKM101 confers Ampicillin resistance).

**Warnings and Precautions:** For Laboratory Use Only

Observe aseptic techniques and established precautions against microbiological hazards throughout all procedures. After use, dispose of all materials according to your institutional biohazard program.

**Storage:** Upon receipt, store plates at 2 – 8°C.

**Procedure:** Streak *E. coli* culture onto each sector of the plate. If streaking from a broth culture with a swab, remove excess broth against the side of the growth vessel prior to streaking.

**Expected Results:**

Sector 1/A: No growth.

Sector 2/B: Growth of all *E. coli* WP2 strains.

Sector 3/C: Growth of pKM101-containing strains only.